

InfraCal 2 Analyzers Measuring Fats, Oils and Grease (FOG) in Wastewater



- Comparable to EPA 1664
- Complies with ASTM D7066
- Battery operable for on-site compliance testing
- Test industrial wastewater prior to release
- Results in under 15 minutes

Industrial and food processing plants that have high levels of fats, oil and grease (FOG) in their effluent face strict wastewater regulations. The InfraCal 2 analyzers provide accurate, easy on-site analysis to help both regulators and industry reduce excessive oil/grease discharges and comply with permit requirements. On-site results are attained in 10-15 minutes, eliminating the wait for remote lab results which can take several

days to a week. The extraction and measurement procedure is simple enough for an operator with minimal training to do the analysis.

The InfraCal 2 is a compact, fixed-filter mid-infrared analyzer with no moving parts. It weighs less than five pounds and can be battery operated. This makes it portable, sturdy and operable in the range of ambient conditions typically found in field environments. The ability to hold multiple calibrations gives pretreatment operators and regulators the ability to measure in the ppm range and the percent range with a single instrument. The infrared technology used in the InfraCal 2 typically correlates well with regulatory methods. The data to the right from an offshore oil rig in the Black Sea compares the InfraCal 2 to the EPA 1664 gravimetric method. The gravimetric method uses 10 times more solvent than the InfraCal 2, takes up to two hours and is not portable. The InfraCal 2 analyzers offer a quick and portable solution for FOG testing.

INFRACAL ANALYZER	NTL-1664
31	33
15	32
17	27
23	29
15	30
20	34
10	12
12	14
12	13
25	24
19	11
13	21
10	21
17	21
17	19
15	19
13	14
13	8

ORDERING INFORMATION

PART NUMBER	ANALYZER MODEL	MINIMUM DETECTION	CALIBRATION OPTIONS	SOLVENT
405-2034	InfraCal 2 ATR-SP	0.3mg/L	Uncalibrated	Hexane, Vertrel MCA, Pentane, Cyclohexane
405-2034-72	InfraCal 2 ATR-SP	0.3mg/L	Factory Cal-Hexane/FOG 6 point set, 10, 25, 50, 100, 150, and 200 mg/L*	Hexane, Vertrel MCA, Pentane, Cyclohexane
405-2035	InfraCal 2 TRANS-SP	0.1mg/L	Uncalibrated	Tetrachloroethylene (perchloroethylene), S-316, Freon
405-2035-40	InfraCal 2 TRANS-SP	0.1mg/L	Factory Cal-S-316/FOG 6 point set-10, 25, 50,100, 150, and 200 mg/L*	S-316

*Note: Lower range calibrations available for oil in water

Spectro Scientific One Executive Drive, Suite 101, Chelmsford, MA 01824-2563 978-431-1120 www.spectrosci.com sales@spectrosci.com An ISO 9001:2008 company Copyright © 2015 Spectro Scientific. All rights reserved. While every effort is made to assure the information in this document is accurate. Spectro Scientific does not accept liability for any errors or mistakes that may arise. Specifications are subject to change without notice. InfraCal 2 FOG_InfraCal2_FOG_2015-10-16